

☒ 2. PL Hi-Plex Ca vs Traditional Low Crosslink Columns For Honey Analysis

A competitor's column, 300x7.8mm

Eluent: Water
Temp: 85°C
Flow Rate: 0.6 ml/min
Column Pressure: 32 bar
Detector: RI

KEY

1. Glucose(4.6K plates/m)
2. Fructose

PL Hi-Plex Ca, 300x7.7mm

Eluent: Water
Temp: 85°C
Flow Rate: 0.6 ml/min
Column Pressure: 15 bar
Detector: RI

KEY

1. Glucose(31.5K plates/m)
2. Fructose

